

# Patch for .NET 3.5 SP1

## Description

The Patch for .NET 3.5 SP1 resolves an issue that causes these software applications to cease functioning when .NET 3.5 SP1 is installed:

- Analyst<sup>®</sup> Reporter software version 1.0
- ChemoView<sup>™</sup> software version 1.0
- Cliquant<sup>®</sup> Food & Beverage Software 1.0
- Cliquant<sup>®</sup> Food & Beverage Software 1.1
- Cliquant<sup>®</sup> Quant Software
- Cliquant<sup>®</sup> Drug Screen & Quant Software
- LightSight<sup>®</sup> software version 2.1
- MRMPilot<sup>™</sup> software version 1.0
- ProteinPilot<sup>™</sup> software version 2.0 and 2.0.1
- AB SCIEX Remote Data Access 2.0 and 2.0.1
- DiscoveryQuant<sup>™</sup> - Analyze software version 2.0.0
- DiscoveryQuant<sup>™</sup> - Optimize software version 2.0.0.1a

## Symptoms

After installing .NET 3.5 SP1, you might experience unexpected crashes or hang-ups while using these software applications.

## Use

The patch automatically finds out which applications are installed on the computer, and replaces a file named rscoree.dll in the applicable installation directories.

## Prerequisites

None

## Installing the Patch for .NET 3.5 SP1

### To install the Patch for .NET 3.5 SP1

- Double-click **Patch.NET3.5SP1.exe** and then follow the instructions in the installation wizard.

### To verify the installation

- Click **Start > Settings > Control Panel > Add/Remove Programs**. The list of currently installed programs should include Patch for .NET 3.5.

**Note:** If you delete the patch by clicking **Change/Remove**, the installation record will be deleted, but rscoree.dll is not deleted from the installation directories.

## Support

Any questions relating to the use of this patch should be directed to Applied Biosystems/MDS Analytical Technologies Technical Support at [ms\\_support@appliedbiosystems.com](mailto:ms_support@appliedbiosystems.com).